## **REAX 2**

**Betriebsanleitung**  
**Instruction Manual**  
**Mode d'Emploi**  
**Instrucciones de**  
**Empleo**  
**Istruzioni per l'uso**

<b>D</b>	DEUTSCH	Seite	3 - 11
<b>E</b>	ENGLISH	page	12 - 20
<b>F</b>	FRANCAISE	page	21 - 29
<b>ES</b>	ESPAÑOL	página	30 - 38
<b>I</b>	ITALIANO	Pagina	39 - 47

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EN 61010

Thank you for purchasing a Heidolph Instruments product. This item has been designed, made and inspected in compliance with DIN EN ISO 61010 for long-term and flawless operation.

**SUMMARY**

**SUMMARY**..... 12  
**STANDARD HARDWARE & OPTIONS**..... 13  
**GENERAL INFORMATION**..... 13  
**SAFETY INFORMATION**..... 14  
**SET-UP**..... 14  
 1. **Set-up**..... 14  
 2. **Electric hook-up**..... 14  
 3. **Secure vessels**..... 14  
   3.1. **Adapter 1**..... 15  
   3.2. **Universal Adapter**..... 15  
   3.3. **Inserting the adapter**..... 16  
**OPERATION AND CONTROLS**..... 16  
**CLEANING & SERVICING**..... 17  
**DISASSEMBLY & STORAGE**..... 17  
**DISPOSAL**..... 18  
**TROUBLESHOOTING**..... 18  
**SPECIFICATIONS**..... 18  
**WARRANTY, LIABILITY & COPYRIGHT**..... 19  
**FAQ / REPAIR WORK**..... 19  
**CE-DECLARATION OF CONFORMITY**..... 20

**Important information**

**Advice about power cord / mains supply**

**Caution: mandatory action**

**Caution: fire- and explosion hazard**

**Advice about maintenance / repair**

## STANDARD HARDWARE & OPTIONS

	Product	quantity	P/N 230/240V 50/60Hz	P/N 115V 50/60Hz
	REAX 2	1	541-21001-00	541-21001-03
	Instruction Manual	1	01-005-001-49	01-005-001-49

### Accessories (optional)

	Item	1st.1.1 P/N
	Adapter for 20 test tubes	549-21000-00
	Universal adapter	549-30000-00

## GENERAL INFORMATION

Unpack your item carefully.  
Inspect for damage and report such damage or missing parts to your supplier right away.

Read your Instruction Manual carefully. Take time to save time when working with your product. Make sure that every user has read and understood the Instruction Manual.

Please store the Instruction Manual in a place easily accessible to every user.

### IF ALL ELSE FAILS, READ THESE INSTRUCTIONS !

A so-called EURO-plug (DIN 49441 CEE 7/VII 10/ 16 A 250 V) is standard on all of the products.  
For the Continental US they feature a US-standard plug (NEMA Pub.No.WDI.1961 ASA C 73.1 . 1961 page 8 15A 125V).

For using the item in a country with deviating outlet / plug systems, we recommend to use approved adapters or to have an electrician replace the standard plug with one suiting your needs.

As shipped, the item features a protective ground wire. When replacing the original plug, make sure to reconnect this protective ground wire in the new plug !

## SAFETY INFORMATION

Please comply with all safety and accident-prevention regulations as in force for laboratory work!

Use extra care when working with flammable substances; refer to safety data sheets.

Use extra care when working in the vicinity of flammable and explosive substances. Motors are of non-sparking type, the item itself however is not explosion-proof.

When connecting your item with your local power supply, please make sure your item is designed for your local voltage; go by the data plate on the item.

Please connect your item with a protective-ground outlet only.

Turn your power switch OFF whenever the item is not used, or before disconnecting the plug.

Repair work is limited to technicians approved by Heidolph Instruments.

Your item needs a solid stand.

Lab bench needs to be of rigid design, and features an anti-skid surface coat.

Before starting the item, make sure all samples are attached safely (must not move while shaking).

## SET-UP

### 1. Set-up

Please locate the shaker on a stable, horizontal surface. For reasons of safety, keep the area around the machine clear of other items.

Be aware of orbital movement of the shaker rotor and vessel set-ups protruding over the standard rotor surface.

### 2. Electric hook-up


Connect mains plug

### 3. Secure vessels

There are two adapters available to meet with a wide range of applications.


### 3.1. Adapter 1

Adapter 1 can hold 20 test tubes with diameters from 10 -18 mm. The vessels are inserted as shown in the illustration.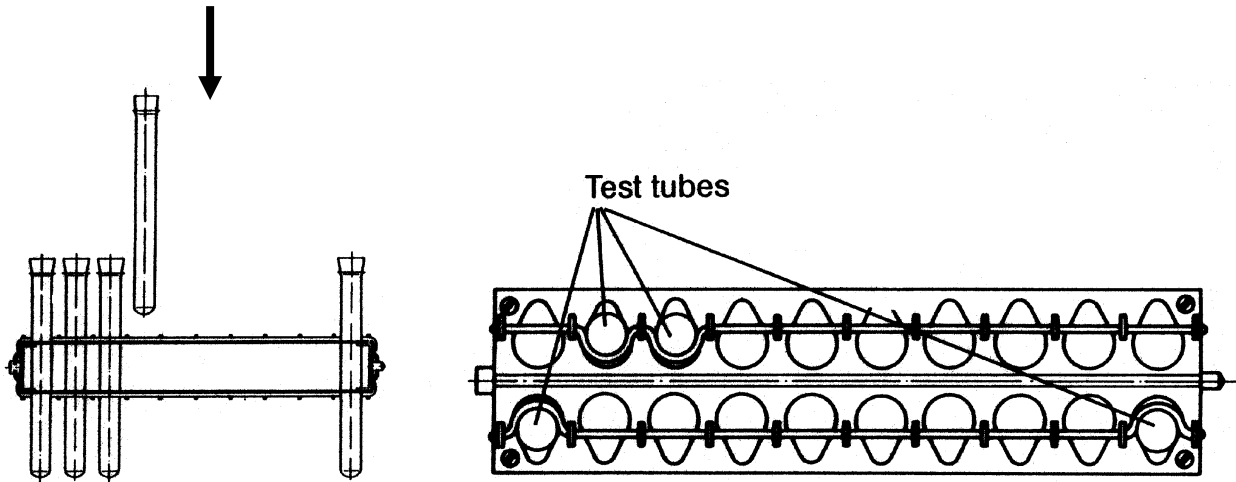


### 3.2. Universal Adapter

The construction of the universal adapter allows to fix different vessels with ground-in stoppers or screwed caps as well as complete test tube racks. To that end, the two quick-action tensioning nuts (H) are loosened by turning them once or twice. When they are in an inclined position, the quick-action tensioning nuts can be moved along the threaded bolt (I). The universal adapter can be opened to a width from 45-160 mm, depending on the case of application. Then the vessels or the test tube rack (K) can be put into the adapter, as shown in the illustration. Next the cover plate (L) is pressed down and secured with the quick-action tensioning nuts (H). (To this end, move the quick-action tensioning nuts along the threaded bolt and tighten them by turning them once or twice.)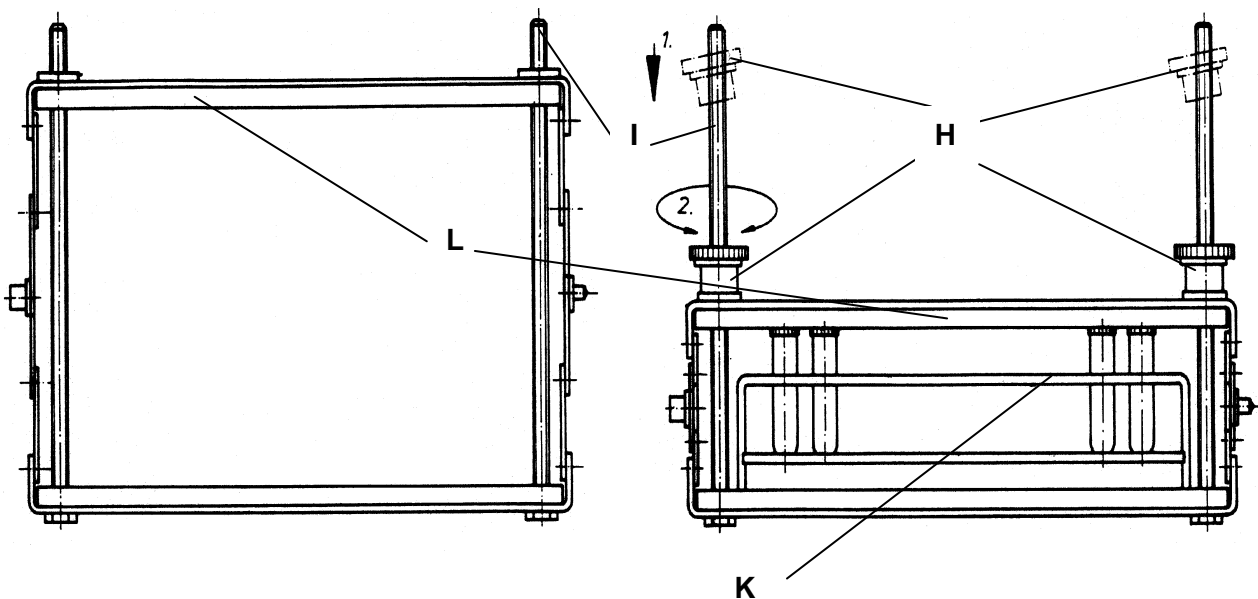

### 3.3. Inserting the Adapter

On one side of the adapter there is a square (C). First this square is inserted into the carrier (D), see illustration. Then the knurled grip (E) is drawn and the pin (F) is placed in the receptacle (G). Finally the knurled grip (E) is brought into its initial position.

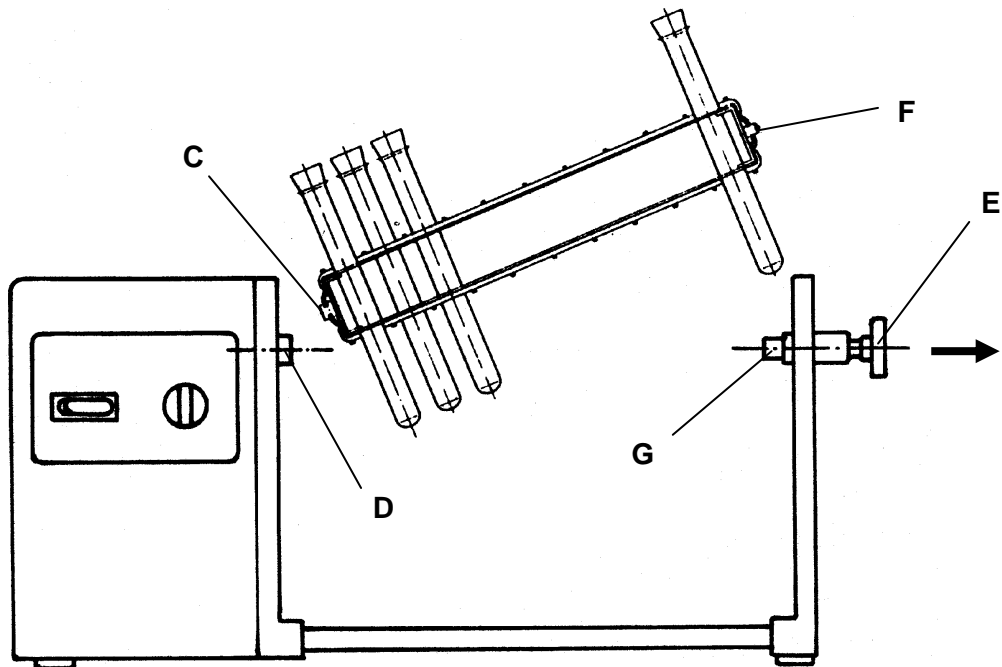## OPERATION AND CONTROLS

Before connecting power cord with mains outlet, make sure that:

- your item is designed for your local voltage and frequency (data plate on item).
- master switch is set to "0" and all controls are in "min." position (turn CCW completely); this way you avoid spilling fluids by too intense shaking action.

Carefully close your vessels and select appropriate shaking intensity (if vessels remain open) to avoid splashes and spillage.


We recommend to start with a low shaking speed and gradually increase frequency to avoid accidental shaking at high speed.

When handling hazardous fluids, make reference to applicable safety information.

- ◆ The appliance is switched on with the main switch (A).
- ◆ Speed Control: The rotation of the adapter is stepless adjustable in the range from 20-100 rpm through the control knob (B). An electronic overspeed monitor provides for constant rotation.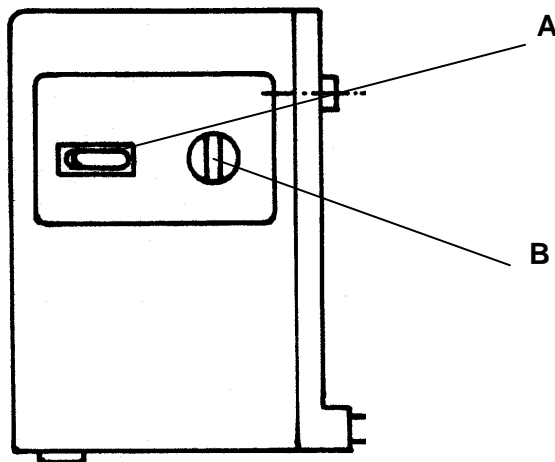

- ◆ The adapter can be turned to the desired position by hand, when the main switch (A) is switched off.

## CLEANING & SERVICING

**Cleaning:** wipe housing clean with a damp cloth (add some sort of mild liquid soap).

### Note

To avoid damage to the surface finish, avoid using chlorine bleach, chlorine-based detergents, abrasive substances, ammonia, rags or cleaning agents containing metal particles.

The item is maintenance-free. Repair work is limited to technicians approved or appointed by HEIDOLPH Instruments. Please call your local HEIDOLPH Instruments Dealer or a HEIDOLPH Instruments Field Representative.

## DISASSEMBLY & STORAGE

### Disassembly

1. Turn item OFF and disconnect mains plug.
2. Remove all of the hardware arranged around the item to ease disassembly.
3. Unload all flasks and beakers from adapter.

### Forward & Store

1. We recommend to store the item and its components in its original box, or a similar container that offers adequate protection against damage in transit. Tape the box securely.
2. Store the item in a dry place.

**Caution**  
Do not jolt or shake the item during transport.

## DISPOSAL

For disposal, please comply with your local or national regulations.  
Split by metal, plastic, etc.  
Packing material to be treated as described above (material split).

## TROUBLESHOOTING

Work on electric and electronic components is limited to qualified personnel.

### Master switch on shaker / mixer won't light

1. check power cord
2. check circuit breakers

### Item won't shake (master switch lighted)

1. Thermal motor circuit breaker triggered by motor overload  
Remedy:  
Wait about 20 minutes, decrease load applied on shaker plate.
2. Mechanical parts broken (humming motor noise) or electronic failure (no motor noise).

## SPECIFICATIONS

### REAX 2

power supply	230/240 V 50/60 Hz / 115 V 50/60 Hz (optional)
input power	27 W
shaker frequency	20 – 100 1/min
mixing movement	overhead rotation
drive	split-pole motor, features electronic speed control; motor overheating protection with automatic reset
adapter	Interchangeable, see accessories
Load-bearing capacity	max. 1 kg, limitations for off-center load
weight	4,8 kg
protection class	IP 21
ambient temperature	0°C to 50°C at 80% rel. humidity Approved for installation in gassing and conditioning cabinets (make reference to temperature limits)
space required	460 mm x 175 mm

## WARRANTY, LIABILITY & COPYRIGHT

### Warranty

Heidolph Instruments warrants that the present product shall be free from defects in material (except wear parts) and workmanship for three years from the date shipped off the manufacturer's warehouse.

Transit damage is excluded from this warranty.

To file for such warranty service, contact Heidolph Instruments (phone:++49-9122-9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty.

Alterations to the present warranty need Heidolph Instruments' consent in writing.

### Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

### Copyright

Copyright in pictures and wording of the present Instruction Manual is held by Heidolph Instruments.

## QUESTIONS / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present Manual, please contact the following address:

For repair services please call Heidolph Instruments (phone: +49 - 9122 - 9920-68) or your local, authorized Heidolph Instruments Dealer.

#### Note

You will receive approval for sending your defective item to the following address:

**Heidolph Instruments GmbH & Co. KG**  
**Lab Equipment Sales**  
**Walpersdorfer Str. 12**  
**D-91126 Schwabach / Germany**  
**Tel.: +49 – 9122 - 9920-74**  
**Fax: +49 – 9122 - 9920-65**  
**E-Mail: Sales@Heidolph.de**

#### Note

If you are based in the United States of America, please contact Heidolph US:

**Heidolph Instruments, LLC**  
**Lab Equipment Sales**  
**2615 River Rd.**  
**Cinnaminson, NJ 08077**  
**Phone: 856-829-6160**  
**Fax: 856-829-7639**  
**E-Mail: heidolph@snip.net**

**Safety Information**

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact *substance*
- take proper protective measure to ensure the safety of our receiving and service personnel
- *mark the pack IAW Hazardous Materials Act*

**CE-DECLARATION OF CONFORMITY**

We herewith declare that the present product complies with the following standards and harmonized documents:

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